

## Smart Clamp™ Cleaning Instructions

Rubber Fab is proud to introduce the next generation of sanitary hinge clamps — the one-piece, completely unitized, Smart Clamp™. With features that include half-turn installation and faster uptime, the benefits of Rubber Fab's Smart Clamp™ far exceed any other sanitary clamp.

**1.** Do not use chlorine bleach based cleaners to clean your Smart Clamp™. We suggest that you use your company's approved cleaner that is not bleach based when cleaning your clamp. This will guarantee a longer product life.

The most important thing to remember when cleaning your Smart Clamp™, either during a CIP, SIP, COP or SOP, is not to get water into the adjustment mechanism of the clamp. Moisture builds up over time and will effect the functionality and life expectancy of the clamp.

Autoclaving is the best option for cleaning the Smart Clamp™. That is completely safe and will keep the Smart Clamp™ in good working order.

In order to keep your Smart Clamp™ in the best working order, there are a couple of things that you should keep in mind when cleaning it.

**2.** Do not submerge your Smart Clamp™ in water or any other liquid. Water or other liquids that have a mineral base could penetrate into the mechanism and leave deposits which could jam the Smart Clamp™.

